


# Vaccine Adjuvants Aluminum Hydroxide Gels (Alum)

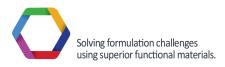



An ABF Ingredients Company

### Vaccine Adjuvants

#### What is an adjuvant?

An adjuvant is a substance that enhances and helps the ability of an antigen to stimulate the immune system. An antigen alone does not stimulate a significant response. Less vaccine antigen (inactive bacteria, virus or toxin) is needed when paired with the right adjuvant.

#### What is VAC?

Aluminum hydroxide gel VAC grades for adsorption (Pseudo-Boehmite form) are used as an adjuvant in human and animal vaccines. It is used to bind negatively charged antigens.

#### Functionality


Its functionality is to adsorb a large range of proteins and inactivated viruses. At pH of preparation of vaccines around 5-7, Pseudo-Boehmite gels have the property to:

- + Adsorb negatively charged antigens
- + Slow the release of an inactivated virus or bacteria
- + Activate and stimulate the immune response
- + Create a more rapid development of high titered and long-lasting antibody responses after primary immunization


#### CUSTOMIZATION

SPI Pharma offers the ability to customize our VAC grades to optimize the performance with your antigen or match your current adjuvant.

Adsorbent properties are related to surface area and surface groups, the hydroxyl groups can accept a proton:

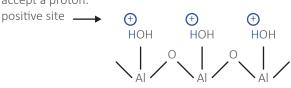

donate a proton:

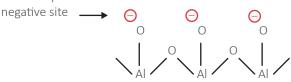


Aluminium Hydroxide Adjuvants surface charge is pH dependant.

Figure 1. Schematic of the structure of the gel. When the surface of the aluminum oxyhydroxide is positive, it attracts negatively charged antigens. Adsorption capacity is determined by surface area and surface charge from the hydroxyl groups.

#### Aluminium Hydroxide gels used as Vaccine Adjuvants:

- + They are called Aluminium Hydroxide but they are not chemically Al(OH)<sub>2</sub>
- + They are poorly crystalline Boehmite (Pseudo-Boehmite): AlO(OH),nH<sub>2</sub>O: Aluminium oxyhydroxide
- + Composed of fibrous primary particles of a few nanometers, Pseudo-Boehmite has adsorption properties and a large surface area (around 500 m2/g)

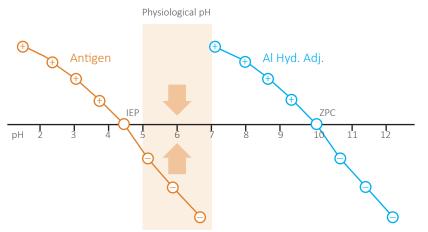

Adjuvants surface charge

Figure 2. At physiological pH around 5.0 – 7.0, Pseudo-Boehmite gels have the property to adsorb negatively charged antigens. Zero point charge (ZPC) is around pH 10.2.

#### EXAMPLE APPLICATIONS

#### Veterinary applications for vaccines for livestock and companion animals (production from inactivated virus)


- + Highly contagious diseases:
  Foot and Mouth Disease, Classical Swine Fever,
  Aujeszky's Disease (Pseudorabies), Infectious Bovine
  Rhinotracheitis (IBR), Avian influenza
- + Swine, Poultry diseases
- + Multiple species diseases:
  Anthrax, Bluetongue, Brucellosis (Brucella abortus),
  Leptospirosis, Paratuberculosis, Q fever, Rabies,
  Rinderpest, West Nile fever

#### SPI'S ADVANTAGES

SPI has two facilities for the production of its VAC products, one in France and the other in US. This allows us to ensure security of supply. Both sites produce under cGMP conditions and are regularly audited by ANSM (French Health Agency), FDA and many customers.

#### Locations where VAC has been sold

- Manufacturing & Sales
- Sales

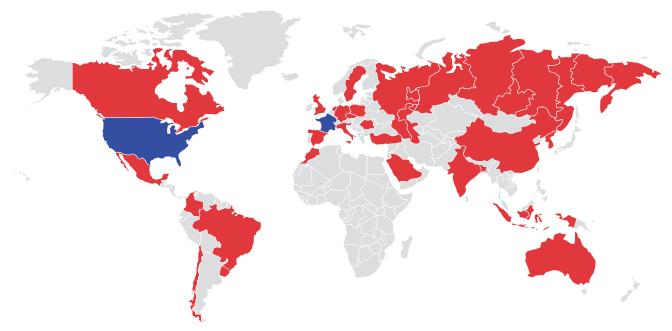



## Vaccine Adjuvants

<b>DESCRIPTION</b> White pumpable suspension, used as a v				
Retest Date: 3 years from the date of ma Storage: Protect from excessive tempera		its original package.		*: around
Grades	Ph. Eur. Specification	VAC 15	VAC 20	VAC 30
Assay Al <sub>2</sub> O <sub>3</sub> % Aluminum stated on the label	90 - 110% Aluminum stated on the label	1.4 - 1.6%	1.9 - 2.1%	2.9 - 3.1%
Corresponding to Al(OH) <sub>3</sub> %		2.14 - 2.45%	2.91 - 3.21%	4.44 - 4.74%
Corresponding to Al %		0.74 – 0.85%	1.00 - 1.11%	1.54 - 1.64%
Protein Binding Capacity, For Info Only		2 mg BSA/mg Al <sub>2</sub> O <sub>3</sub> *		
Adsorption Power (Ph. Eur. Test)	Conforms No detectable BSA in solution	Conforms No detectable BSA in solution	Conforms No detectable BSA in solution	Conforms No detectable BSA in solution
Point of Zero Charge, For Info Only		10.2	10.2	10.2
Sedimentation (Ph. Eur. Test) @ 5mg Al/mL	NMT 5 mL	NMT 5 mL	NMT 5 mL	NMT 5 mL
Viscosity, For Info Only		NMT 500 cps (Brookfield V3/30)	NMT 1000 cps (Brookfield V3/30)	NMT 1500 cps (Brookfield V3/30)
Nitrate (NO <sub>3</sub> ) (Ph. Eur. Test)	NMT 100 ppm	NMT 100 ppm	NMT 100 ppm	NMT 100 ppm
Ammonium (NH <sub>4</sub> ) (Ph. Eur. Test)	NMT 50 ppm	NMT 50 ppm	NMT 50 ppm	NMT 50 ppm
Sulfate (Total SO <sub>4</sub> ) (Ph. Eur. Test)	NMT 0.5%	NMT 0.5%	NMT 0.5%	NMT 0.5%
Chloride (Cl) (Ph. Eur. Test)	NMT 0.33%	NMT 0.33%	NMT 0.33%	NMT 0.33%
Heavy Metals (as Pb) (Ph. Eur. Test)	NMT 20 ppm	NMT 20 ppm	NMT 20 ppm	NMT 20 ppm
Arsenic (As) (Ph. Eur. Test)	NMT 1 ppm	NMT 1 ppm	NMT 1 ppm	NMT 1 ppm
Iron (Fe) (Ph. Eur. Test)	NMT 15 ppm	NMT 15 ppm	NMT 15 ppm	NMT 15 ppm
pH (Ph. Eur. Test)	5.5 – 8.5	5.5 - 8.5	5.5 – 8.5	5.5 – 8.5
Endotoxins (Ph. Eur. Test)	LT 5 IU	LT 5 IU	LT 5 IU	LT 5 IU
Standard Packaging		50kg plastic drum	25kg Jerrycan	50kg plastic drum
Sterilization		All products can be sterilized.		

Americas SPI Pharma, Inc. Rockwood Office Park 503 Carr Rd., Suite 210 Wilmington, DE 19809

Contact Us: salesinfo@spipharma.com Europe/Middle East/ Africa SPI Pharma SAS Chemin du Vallon du Maire 13240 Septemes-Les Vallons France Asia/Pacific SPI Pharma, Inc – India Branch 21 B, Veerasandra Industrial Area Hosur Road, Bangalore – 560100 Karnataka, India Australia Distribution Company Anzchem 1 Braidwood Street Enfield NSW 2136 Australia www.spipharma.com



© SPI Pharma 2017. All trademarks are the property of SPI Pharma. The information contained in this document is proprietary to SPI Pharma and may not be used or disseminated inappropriately. The information and recommendations contained herein are to the best of SPI Pharma, Inc.'s knowledge reliable and accurate. Any recommendations are made without warranty, either implied or expressed, due to the variations in equipment, conditions, and methods which may be used in commercially processing the products. No warranties of any kind are made, express or implied, including those of merchantability and fitness for particular purpose, other than the products conform to current standard specifications. SPI Pharma, Inc. makes no warranty that the use of the products or formulations provided by SPI Pharma, Inc. will not infringe any trademark, trade name, copyright, patent or other rights held by any third party when used in customer's application. SPI Pharma, Inc. shall not be liable for loss of profit or for incidental, special or consequential loss or damages.

Order No. SPI-VAC-VAC-0100-10201701 10-2017 | All rights reserved